
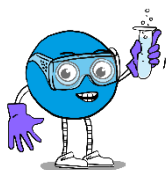


Peptide Solubilization Protocol

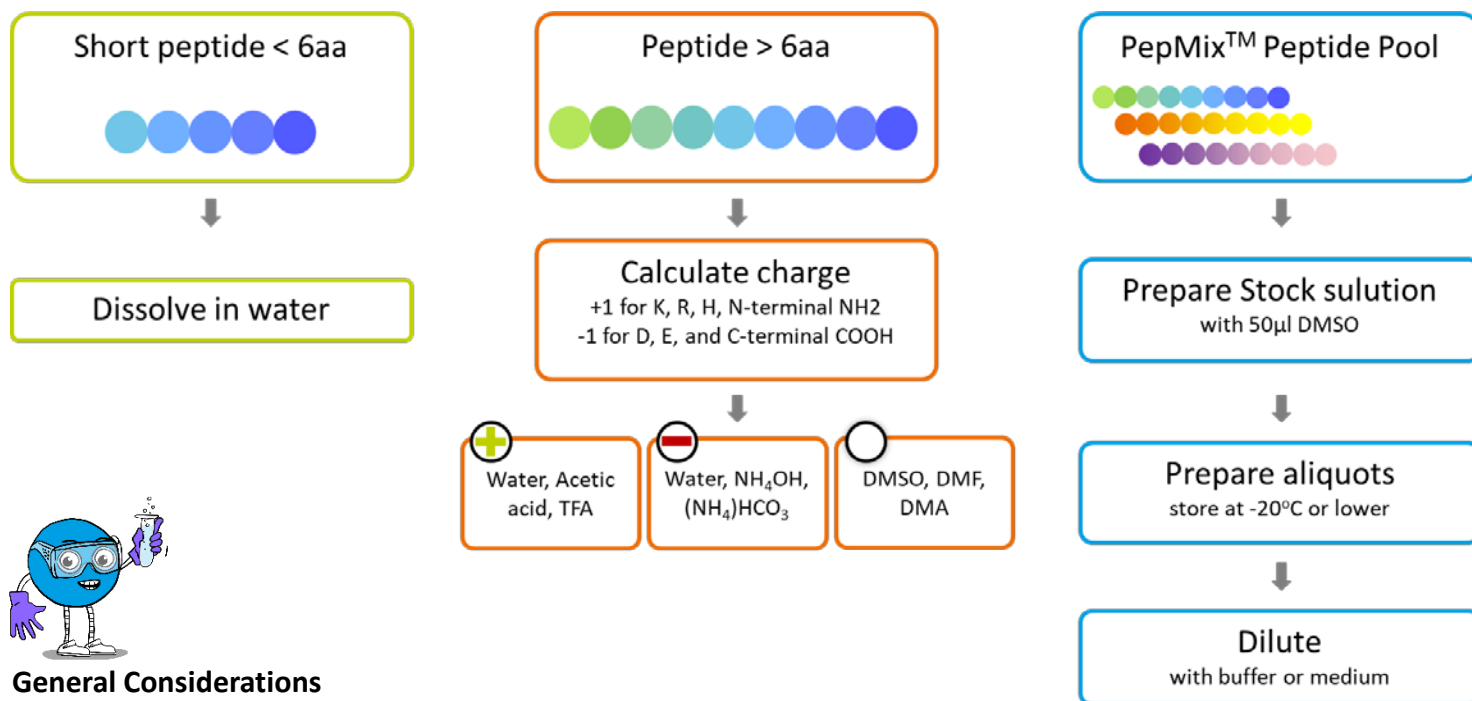

General Considerations

- Always test solubility first with a tiny amount of peptide.
- Do not to exceed 1% (v/v) DMSO in your final solution if you are using your peptides in cellular assays.
- DMSO should not be used with peptides containing methionine or cysteine as it may oxidize the side chain. Use DMF instead.
- Vortex or sonicate after each step.
- Centrifuge your peptide solution before use to precipitate any undissolved peptide residues.

Visit the
Peptide Expert!

[https://www.jpt.com/
peptide@jpt.com](https://www.jpt.com/peptide@jpt.com)

**We produce over 1
million peptides per year!**

ISO 9001:2015 certified

Peptide
Synthesis
Service



Peptide
Solubility
Testing

